



# FUNGAL ASSESSMENT BY CULTURE AND RT-PCR ASSAY IN A WASTE SORTING PLANT



## CORIOLIS AIR SAMPLERS USED TO PREVENT WORKERS EXPOSURE TO POLLUTANTS

Polytechnic Institute of Lisbon, Lisbon, Portugal

### CONTEXT

In the management of **solid waste**, **pollutants** over a wide range are released with different routes of exposure for workers. The potential for synergism among the pollutants raises concerns about potential adverse **health effects**, and there are still many uncertainties involved in exposure assessment.

In this study [1], **molecular real-time polymerase chain reaction (RT-PCR)** was used to assess the presence of three potential **pathogenic/toxigenic** fungal species: *Aspergillus flavus*, *A. fumigatus*, and *Stachybotrys chartarum* from air samples collected with a **Coriolis  $\mu$**  in a waste-sorting plant. These results were compared with the cultivable fungal air contamination obtained by conventional method (culture-based).

### MATERIALS

- **Coriolis  $\mu$** , sterile cones, 10mL sterile phosphate-buffered saline with 0.005% TritonX100.
- **Conventional method:** Millipore air Tester (Millipore), sterile petri dishes, malt extract agar (MEA) with chloramphenicol (0.05%).





# FUNGAL ASSESSMENT BY CULTURE AND RT-PCR ASSAY IN A WASTE SORTING PLANT

## PROTOCOL

- Sampling at five different workstations inside the facilities and also outdoors as a reference.
- Coriolis air sampling: 300L/min, 1min. Liquid air sample analyzed by RT-PCR (Specific primers and TaqMan probes were used to detect the presence of each specie by Real Time PCR in iQ Real Time Detection System (Bio-Rad)).
- Culture-based analysis (Conventional method): 50L collected by impaction method (140 L/min). Incubation of agar plates for 5 to 7 days at +27 °C.

## RESULTS

Penicillium sp. were the most common species found at all plant locations by conventional methodology. Pathogenic/toxigenic species (*A. fumigatus* and *S. chartarum*) were detected at two different workstations by RT-PCR but not by culture-based techniques (Table 1).

Our results illustrated the advantage of combining both conventional and molecular methodologies in fungal exposure assessment.

Together with microbial volatile organic compounds (MVOC) analyses in indoor air, data obtained allow for a more precise evaluation of potential health risks associated with bioaerosol exposure. Consequently, with this knowledge, strategies may be developed for effective protection of the workers.

Samples	Fungi quantification (CFU/m <sup>3</sup> )	Pathogenic/toxigenic species detection*	
		Culture	Real-time PCR
LDPE 1	<i>Penicillium</i> (full plate)	Not detected	
LDPE 2	<i>Penicillium</i> (full plate)	Not detected	
Paper 1	<i>Penicillium</i> (full plate)	Not detected	<i>A. fumigatus</i>
Paper 2	<i>Penicillium</i> (full plate)	Not detected	
Hall	<i>Penicillium</i> (full plate), 40 <i>Rhizopus</i>	Not detected	<i>A. fumigatus</i> <i>S. chartarum</i>
Outdoor	320 <i>Penicillium</i> , 20 <i>A. clavatus</i> , 20 <i>Chrysosporium</i>	Not detected	

\* *A. flavus*, *A. fumigatus*, *S. chartarum*

[1] J. Malta-Vacas et al, Fungal and Microbial Volatile Organic Compounds Exposure Assessment in a Waste Sorting Plant, Journal of Toxicology and Environmental Health, Part A: Current Issues Volume 75, Issue 22-23, 2012, doi:10.1080/15287394.2012.721175




# FUNGAL ASSESSMENT BY CULTURE AND RT-PCR ASSAY IN A WASTE SORTING PLANT

## CONCLUSION

- Molecular methods are essential for the detection of pathogenic/toxigenic species, especially in highly contaminated environments, where their presence can be masked by other predominant species.
- Coriolis  $\mu$  air sampler offers the advantage of combining both conventional (culture-based) and molecular analysis (RT-PCR) to assess fungal air contamination in waste sorting plant.



## DISCOVER OUR COMPREHENSIVE RANGE OF SOLUTIONS

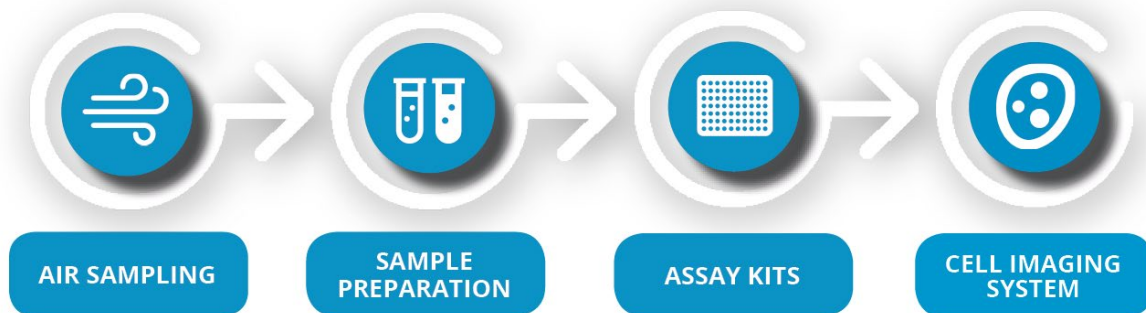

**Join the Bertin Instruments community!**

1500+ documents available online on the Application center

[www.bertin-instruments.com/application-center/](http://www.bertin-instruments.com/application-center/)